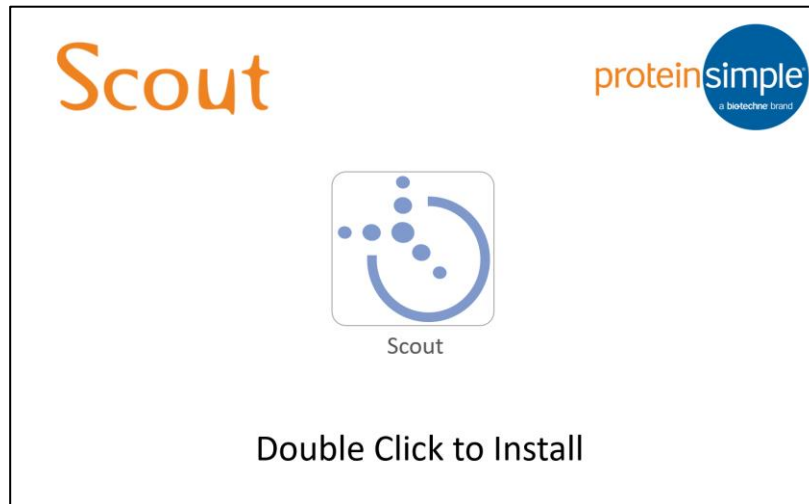


**2.1.10**

The 2.1.10 release is a signed macOS application. Install by opening the dmg file and then double-clicking on the installation application. The contents of the dmg will look like:

**2.0.5.beta**

If you receive a security warning when attempting to install the 2.0.5.beta macOS version of Scout:

1. In the Finder, locate the app you want to open.

Don't use Launchpad to do this. Launchpad doesn't allow you to access the shortcut menu.

2. Press the Control key and click the app icon, then choose Open from the shortcut menu.
3. Click Open.

The app is saved as an exception to your security settings, and you can open it in the future by double-clicking it just as you can any registered app.

Note: Another way to grant an exception for a blocked app is to click the "Open Anyway" button on the Security pane of System Preferences. This button is available for about an hour after you try to open the app.

To open this pane, choose Apple menu > System Preferences, then click Security & Privacy.

## Scout macOS installation notes

For installation on macOS Sierra (10.12), please copy the .mac64 installation program from inside of the downloaded .dmg file to your desktop (or other disk location) before running the installation program. The installation may fail if you attempt to run the .mac64 program directly from inside the downloaded .dmg file. See the figure below.

